

Glucose Glutamic Acid Single-Use Vial For BOD testing:

The instructions are simple... Shake thoroughly, remove the celloseal covering the cap, unscrew the cap, pour contents, and rinse emptied vial at least 2 times with dilution water into the check standard solution. Each vial contains 6mL of the APHA specified concentration of Glucose Glutamic Acid (150mg/L of Glucose, 150mg/L of Glutamic Acid).

We know you will find that this single-use vial improves test accuracy, reduces cost, and eliminates the risk and waste of GGA contamination.

If you have any questions concerning use or formulation - or if you run into any difficulties – please contact one of our Chemists at:

LabChem Inc

Tel: 412-826-5230 Email: mattp@labchem.com

> Catalog Number LC14845-246 is the 24 vial pack Catalog Number LC14845-486 is the 48 vial pack

Enjoy!